IMMB-06-03-05-22-EN-PL

Start up Guide

Szybki start

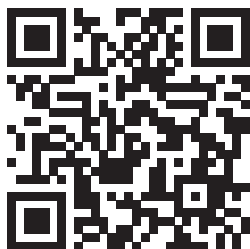
Precision Balances

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- WLC A1
- WLC A2
- WLC A1/C/2
- WLC A2/C/2

To view a complete User Manual go to the website or scan QR code:

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
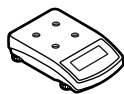
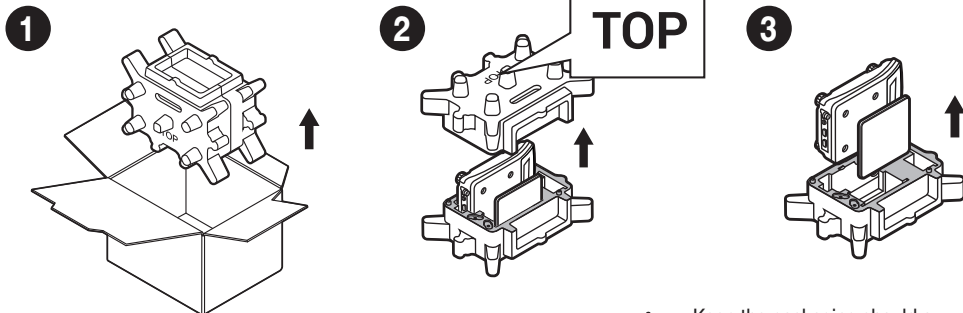
3. RECOMMENDATIONS AND WARNINGS

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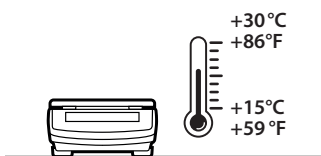
Operate the device in a room where the temperature ranges between 15–30 °C (59–86 °F) and where the relative humidity is below 80%. Any changes in temperature should be gradual and very slow.


Place the balance on a solid surface to ensure stability. To obtain stable and repeatable weighing results, an anti-vibration table is recommended.

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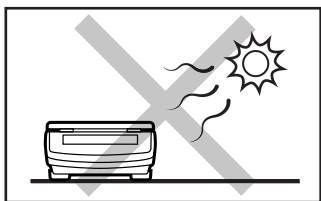
Place the balance away from heat sources. Avoid exposing the balance to the sunlight.

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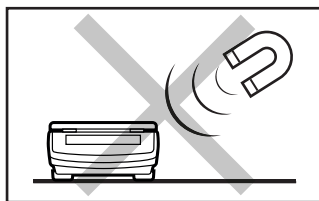
Avoid exposing the balance to a magnetic field. Do not weigh magnetic substances.

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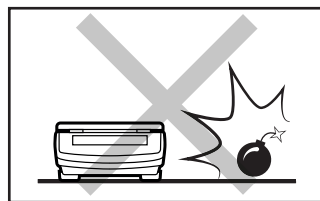
Do not place the balance in a hazardous area. Do not weigh explosive materials.

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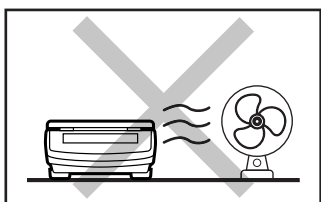
Avoid air drafts and air movements at the workstation.

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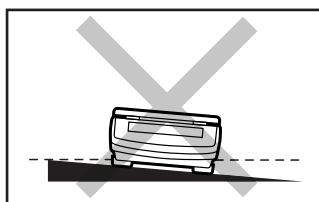Make sure that the balance is placed on an even surface.

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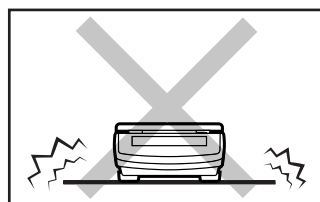Do not place the balance on an unstable ground exposed to shocks and vibrations.

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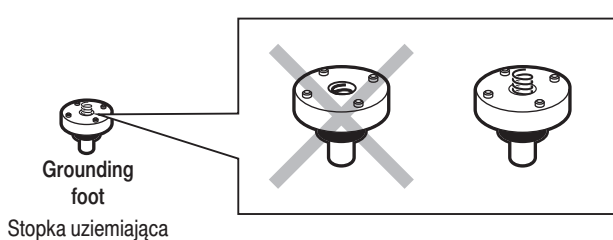
4. ACTIVITIES TO BE DONE PRIOR TO OPERATION

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4.1. Grounding spring check

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
Check whether the grounding spring takes its intended location. Make sure that the grounding spring juts slightly out of the hole.

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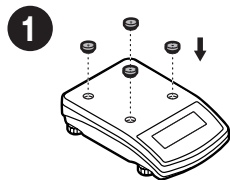
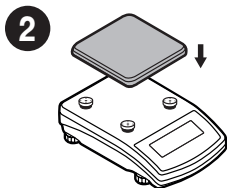
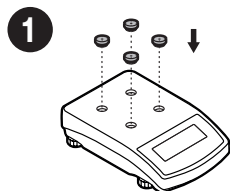
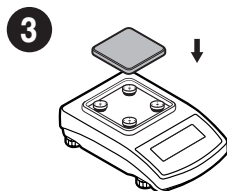
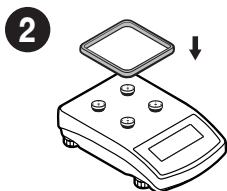
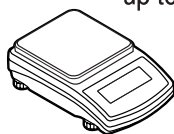
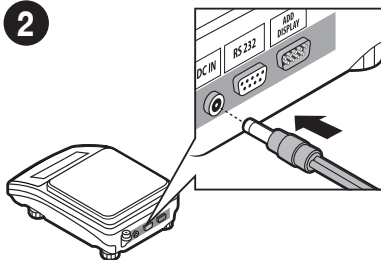
2

3

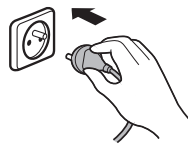Prior to switching the device on, let it reach room temperature (estimated stabilisation time: up to 8 hours).

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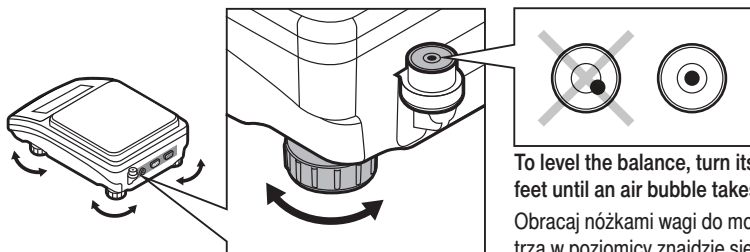
6.2. Levelling

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
To level the balance, turn its feet and keep turning the feet until an air bubble takes the central position.

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6.3. Switching the balance on

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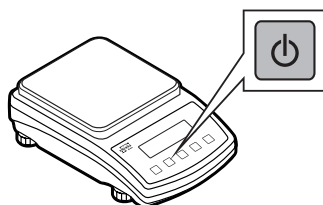To switch the device on, press ON / OFF key.

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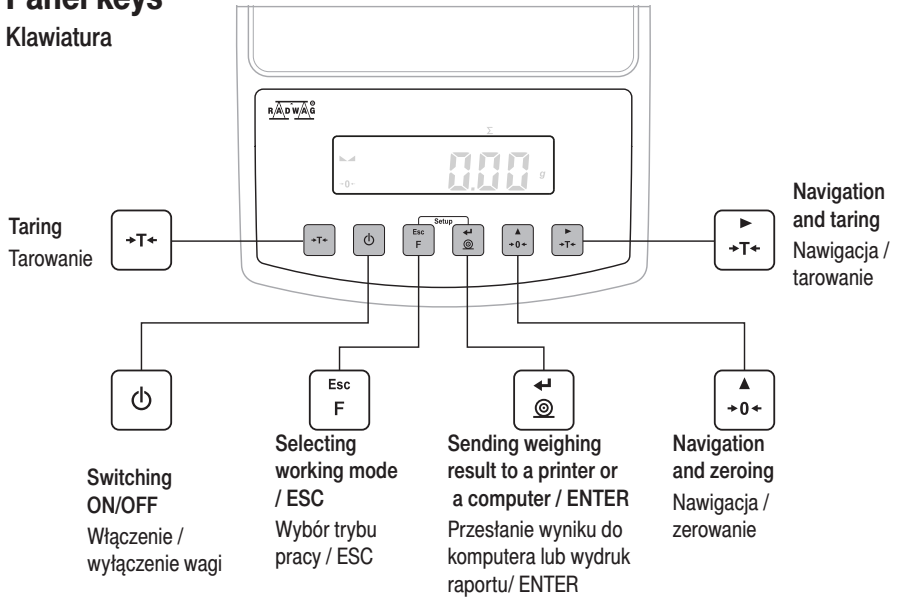
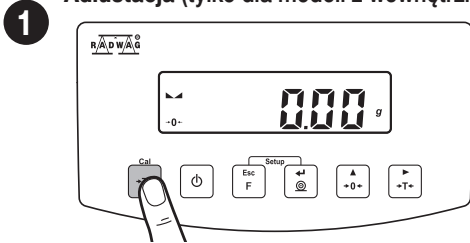
8. BASIC OPERATIONS

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
8.1. Adjustment (for models with an internal adjustment only)

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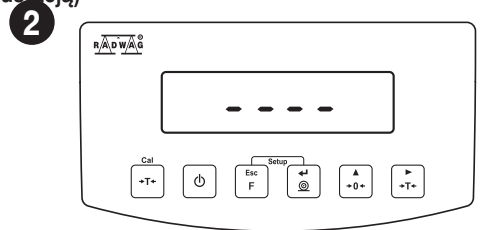8.2. Adjustment (for models with an external adjustment only)

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External adjustment is carried out using an external mass standard of the right accuracy and weight value, which value depends on balance type and capacity.

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
The adjustment mass standard does not come with the balance. Prior to adjustment prepare the right mass standard. You are recommended to use a mass standard of class F1 or F2. View the table and select the mass standard you need.

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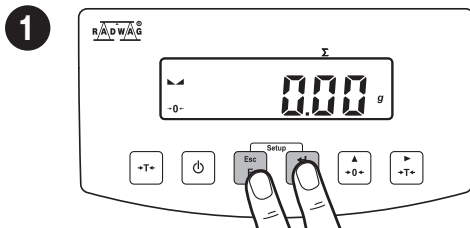
Simultaneously press **Esc F** + **Setup** keys to enter home screen.

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
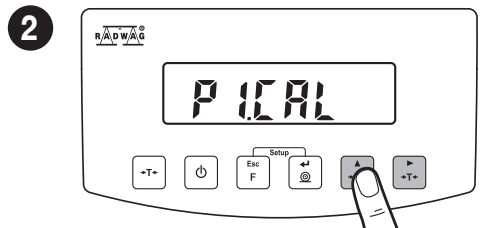
* - N/A in the case of a non-verified instrument

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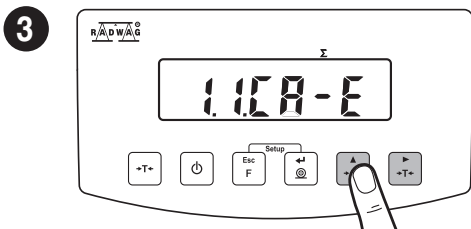
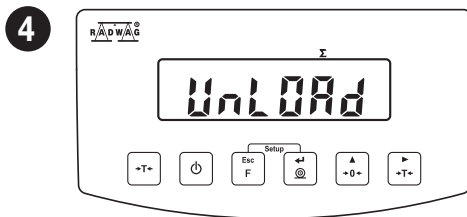
Wait for completion of the start mass determination process.

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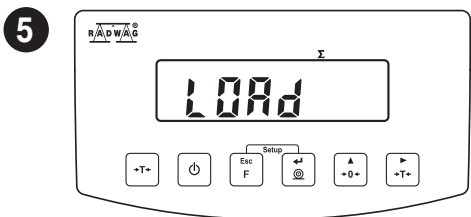
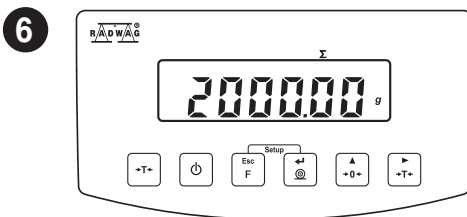
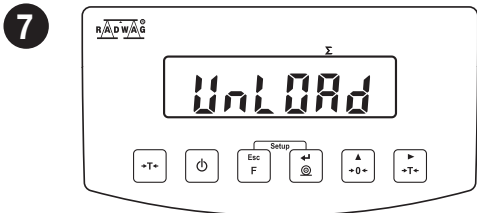
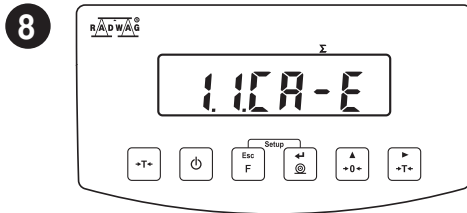
8.3. Zeroing

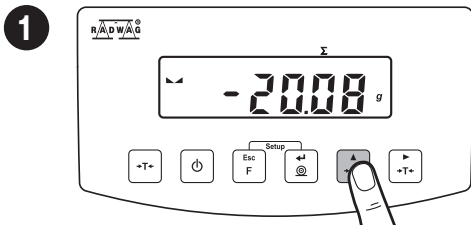Zerowanie

Make sure that the weighing pan is empty and press **+0-** key.

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
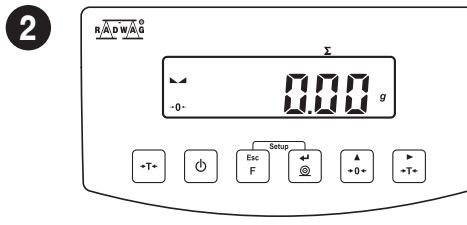
8.4. Taring

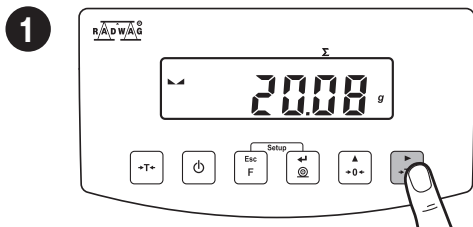Tarowanie

With loaded weighing pan: upon stabilisation of the weighing result, press **-T-** key.

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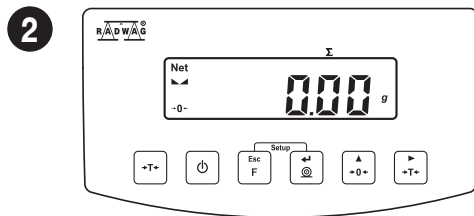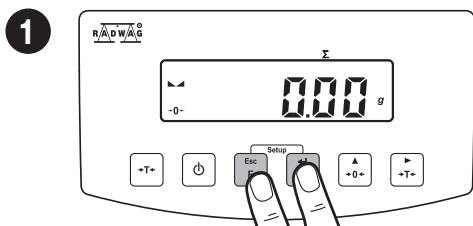
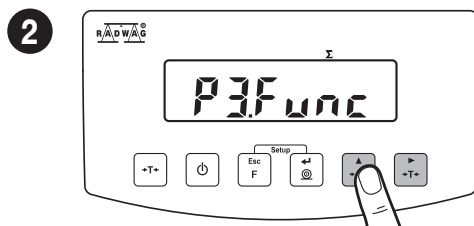
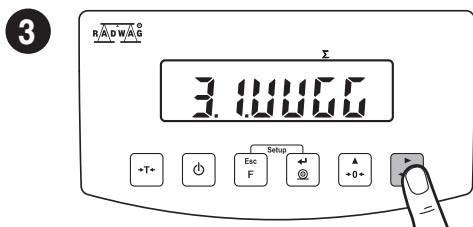
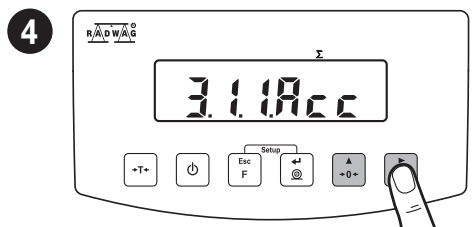
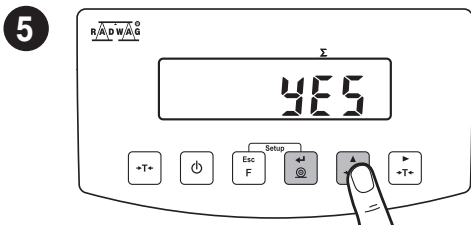
Press key to select “YES” value, next press key in order to save settings.

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To configure other working modes, follow the above procedure.

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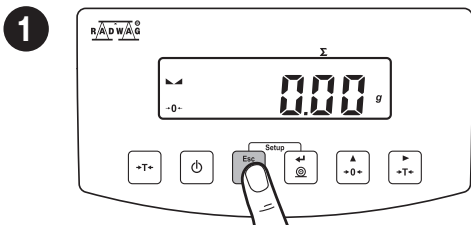
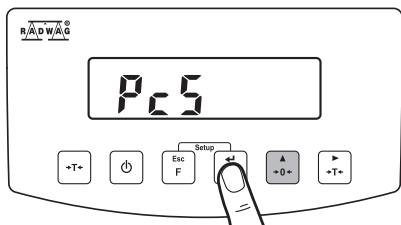
2

Using key select working mode and press .

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The menu features the following working modes:

- Parts counting
- Percent weighing
- Checkweighing
- Animal weighing
- Totalising
- Peak hold
- Weighing

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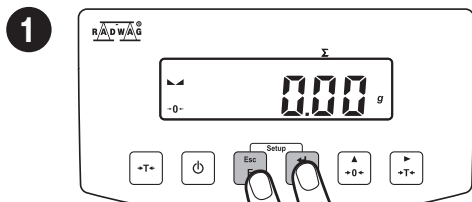
Detailed overview of the settings of the working modes is to be found in balance's main user manual.

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
8.7. Weighing unit selection

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
Simultaneously press + keys to enter home screen.

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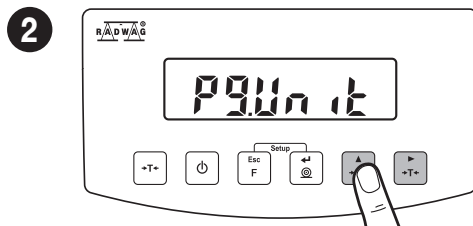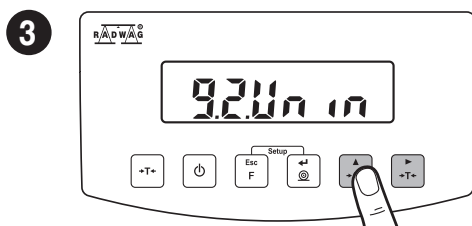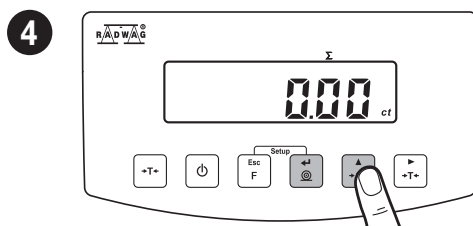
Using key select weighing unit and press key to confirm.

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The balance enables selecting the following weighing units:

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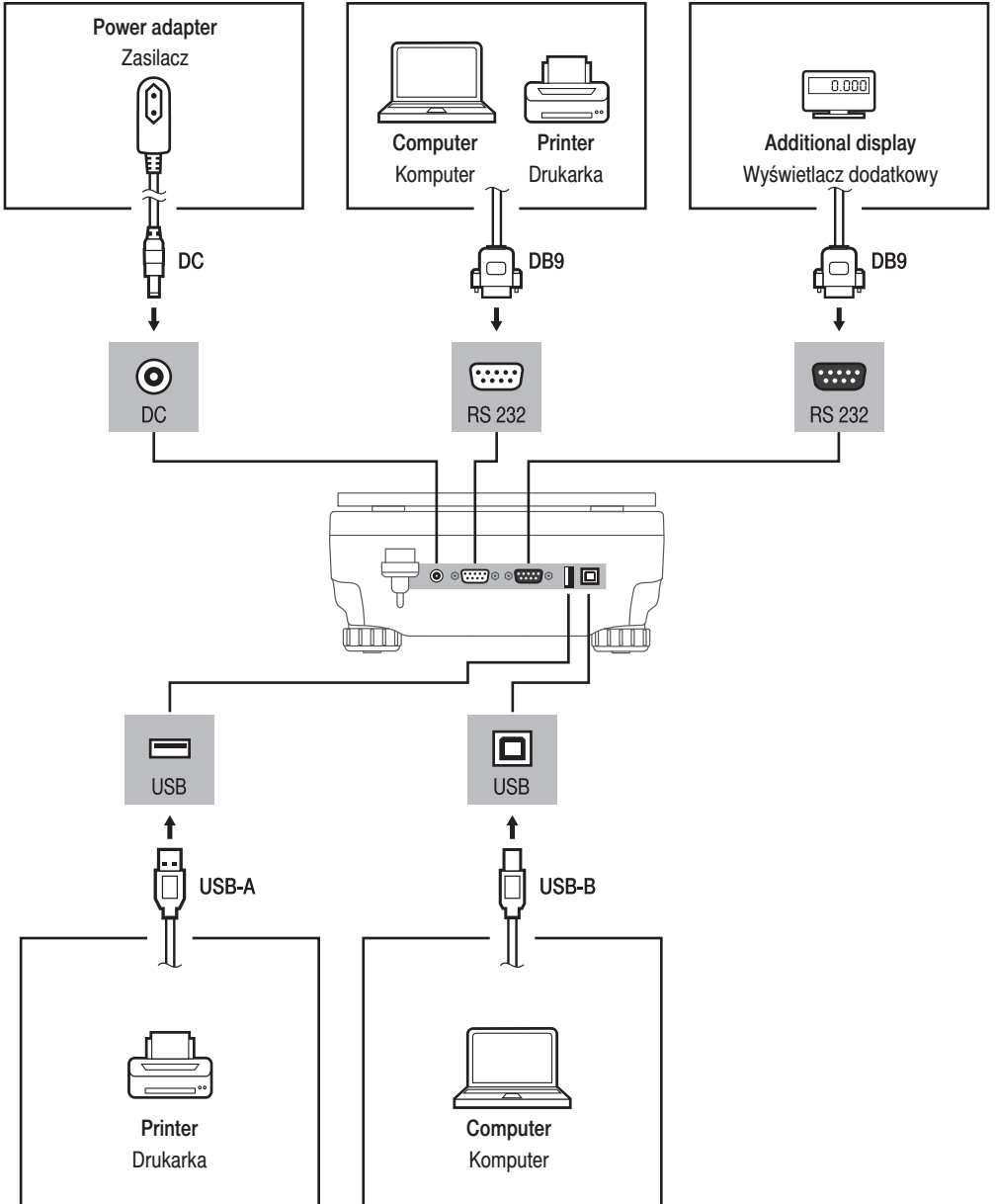
- g (gram)
- kg (kilogram)
- ct (karat)
- lb (funt)*

*(lb) pound unit unavailable for verified balances.

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
10. DEVICE CLEANING

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
10.1. General recommendations

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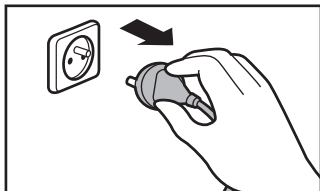
Prior to cleaning, it is necessary to disconnect the balance from the mains.

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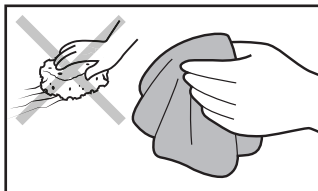Use soft cloths made of microfiber, natural fiber or man-made fiber. Avoid using abrasive cloths or cloths that might scratch the surface!

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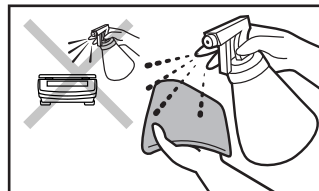Apply the cleanser onto the cloth first. Avoid applying the cleanser directly onto the device!

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Do not use aggressive cleaning agents (e.g. solvents, chlorine preparations, corrosive chemicals, bleach).

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Neither use substances of sharp structure nor intended for scrubbing.

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Make sure that neither dust nor liquid gets inside the weighing system (device inside).

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- **GLASS PANES AND COMPONENTS** can be cleaned using window cleaner.
- **WEIGHING PANS AND STAINLESS STEEL OR ALUMINIUM COMPONENTS** can be cleaned using a cloth slightly soaked in a solution of water and gentle cleanser (e.g. soap or dishwashing liquid).
- **OPERATION PANEL AND HOUSING** can be cleaned using a cloth slightly soaked in a solution of water and gentle cleanser (e.g. soap or dishwashing liquid).

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