IMMB-08-03-06-24-EN-PL

Start up Guide

Szybki start

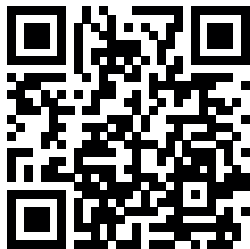
Precision balances

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- WTC

To view a complete User Manual go to the website or scan QR code:

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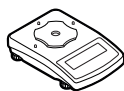
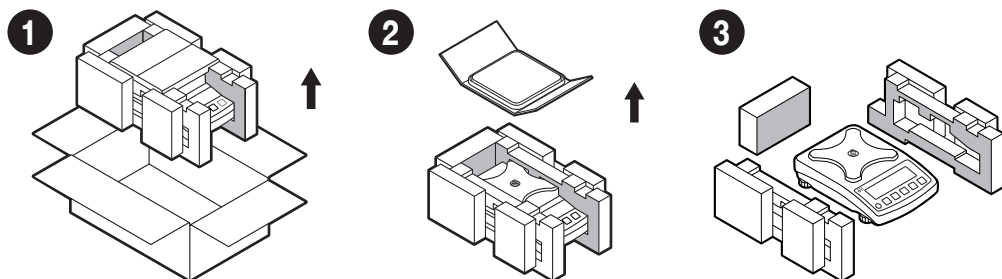
2. UNPACKING

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
Keep the packaging should a warranty claim or service be required.

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
3. RECOMMENDATIONS AND WARNINGS

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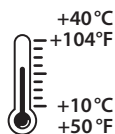
Operate the device in a room where the temperature ranges between 10 – 40 °C (50–104 °F) and where the relative humidity is below 80%. Place the balance on a solid surface to ensure stability. To obtain stable and repeatable weighing results, an anti-vibration table is recommended.

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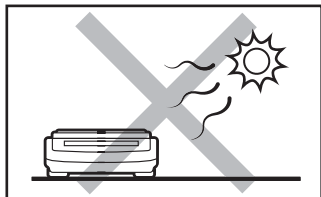
Place the balance away from heat sources. Avoid exposing the balance to the sunlight.

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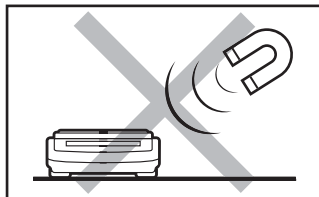
Avoid exposing the balance to a magnetic field. Do not weigh magnetic substances.

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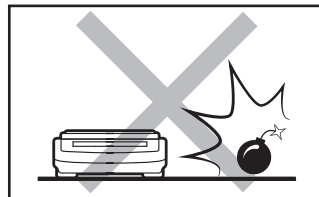
Do not place the balance in a hazardous area. Do not weigh explosive materials.

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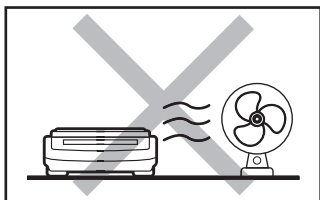
Avoid air drafts and air movements at the workstation.

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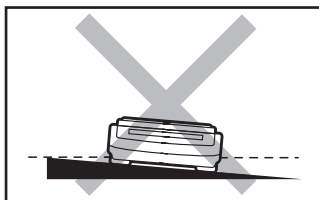Make sure that the balance is placed on an even surface.

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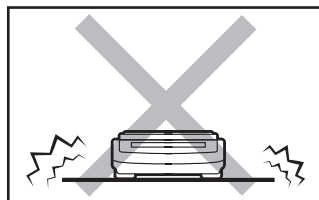Do not place the balance on an unstable ground exposed to shocks and vibrations.

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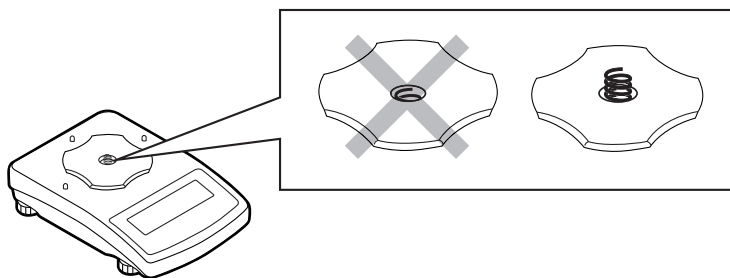
4. ACTIVITIES TO BE DONE PRIOR TO OPERATION

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4.1. Grounding spring check

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

Check whether the grounding spring takes its intended location. Make sure that the grounding spring juts slightly out of the hole.

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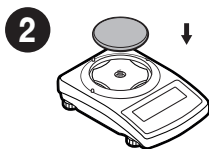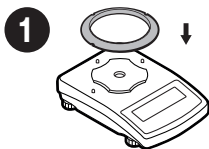
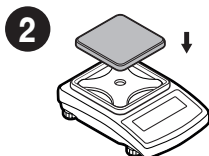
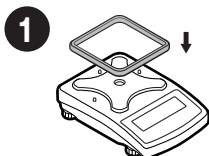
6. PREPARING FOR WORK

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6.1. Temperature stabilization

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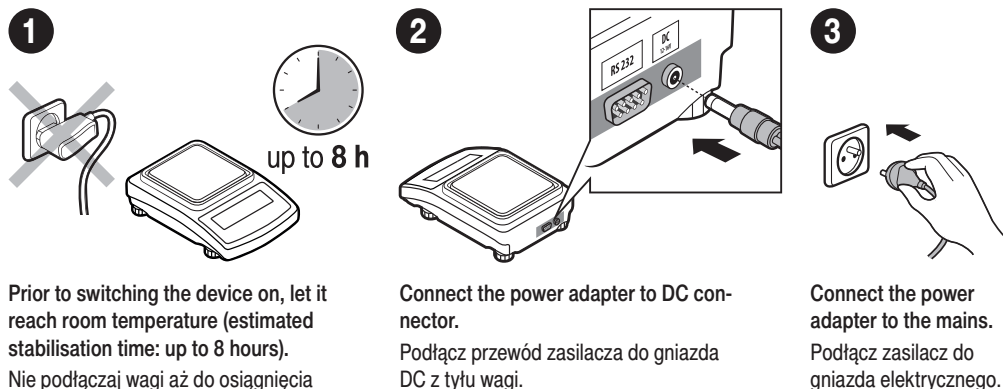Prior to switching the device on, let it reach room temperature (estimated stabilisation time: up to 8 hours).

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Connect the power adapter to DC connector.

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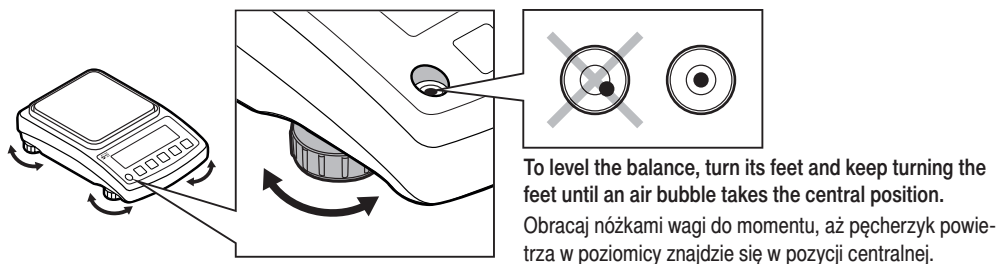
Connect the power adapter to the mains.

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
6.2. Levelling

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6.3. Switching the balance on

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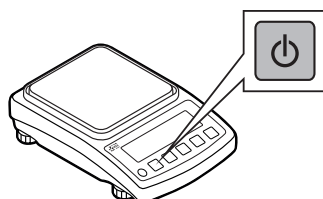To switch the device on, press ON / OFF key.

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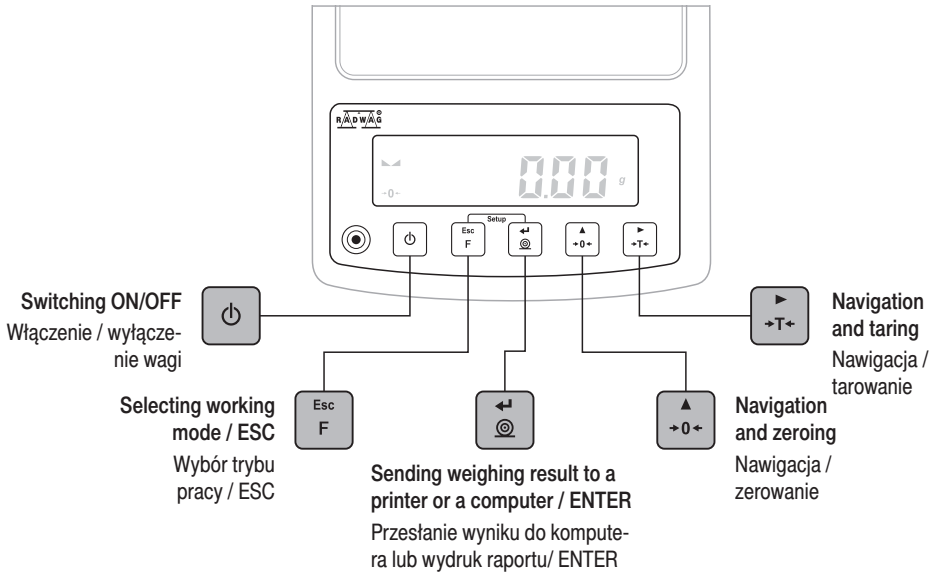
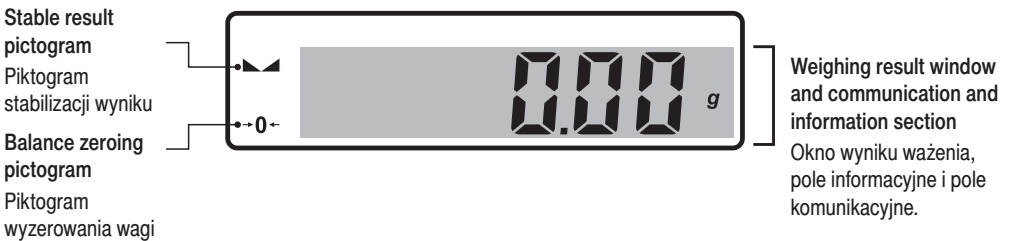
8.1. Adjustment

Adiustacja

External adjustment is carried out using an external mass standard of the right accuracy and weight value, which value depends on balance type and capacity.

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
The adjustment mass standard does not come with the balance. Prior to adjustment prepare the right mass standard. You are recommended to use a mass standard of class F1 or F2. View the table and select the mass standard you need.

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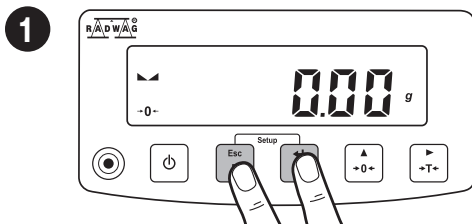Simultaneously press + keys to enter home screen.

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

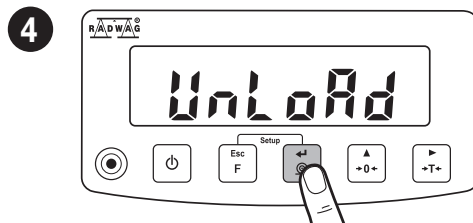
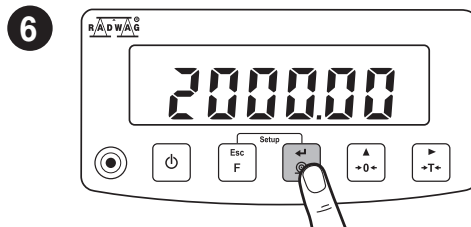
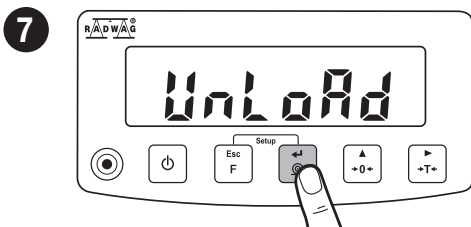
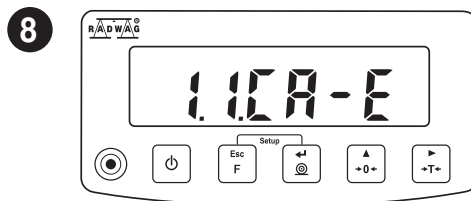
8.3. Taring

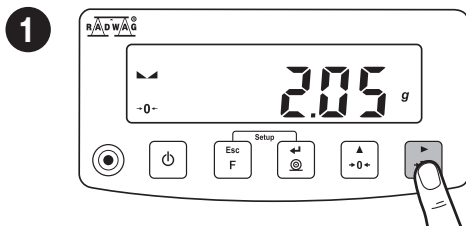Tarowanie

With loaded weighing pan, upon result stabilization, press **-T-** key.

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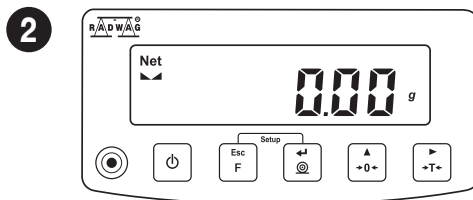
The balance has been tared.

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
8.4. Working modes accessibility

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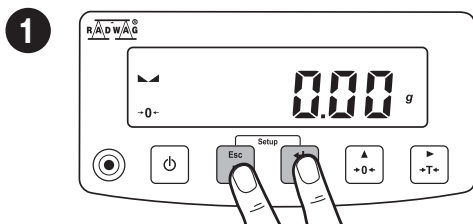
Simultaneously press **Esc F** + **Setup** keys to enter home screen.

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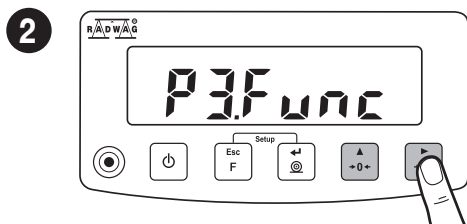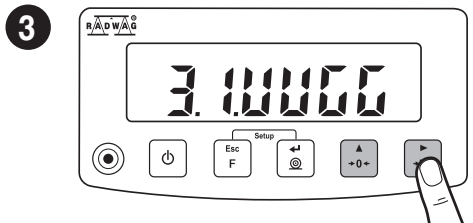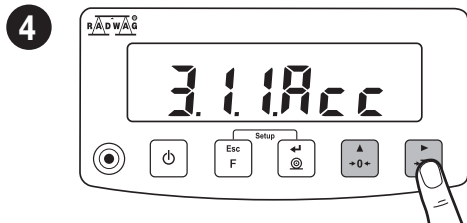
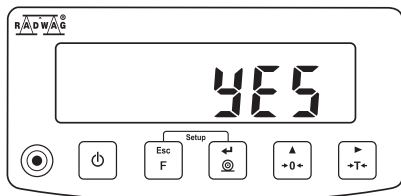
5

Using key select "YES" value and press key to confirm.

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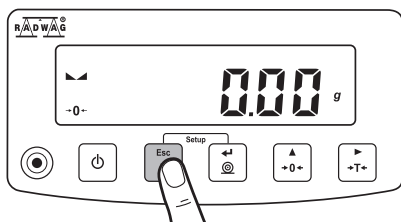
To configure other working modes, follow the above procedure.

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
Using key enter working modes menu.

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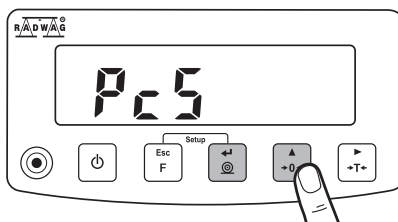2

Using key select working mode and press .

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The menu features the following working modes:

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Detailed overview of the settings of the working modes is to be found in balance's main user manual.

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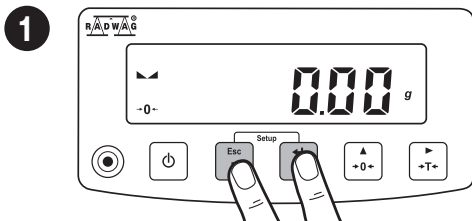
8.6. Weighing unit selection

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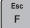
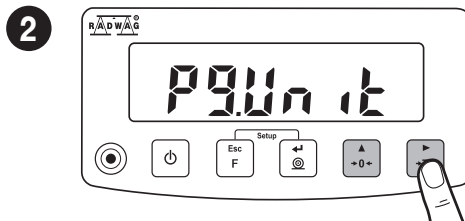
When the main unit is g (gram), the user has the following units: g (gram), ct (karat), lb (pound).*

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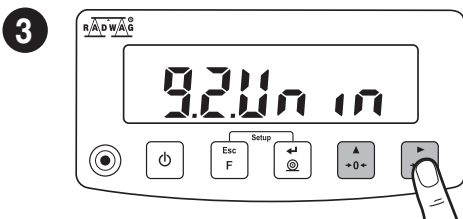
1 Simultaneously press   keys to enter home screen.

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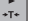
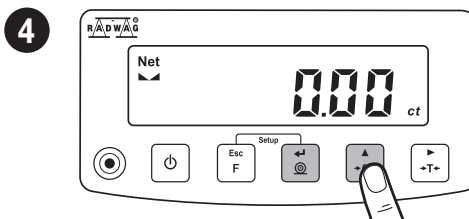
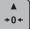
4 Using  key select weighing unit and press  key to confirm.

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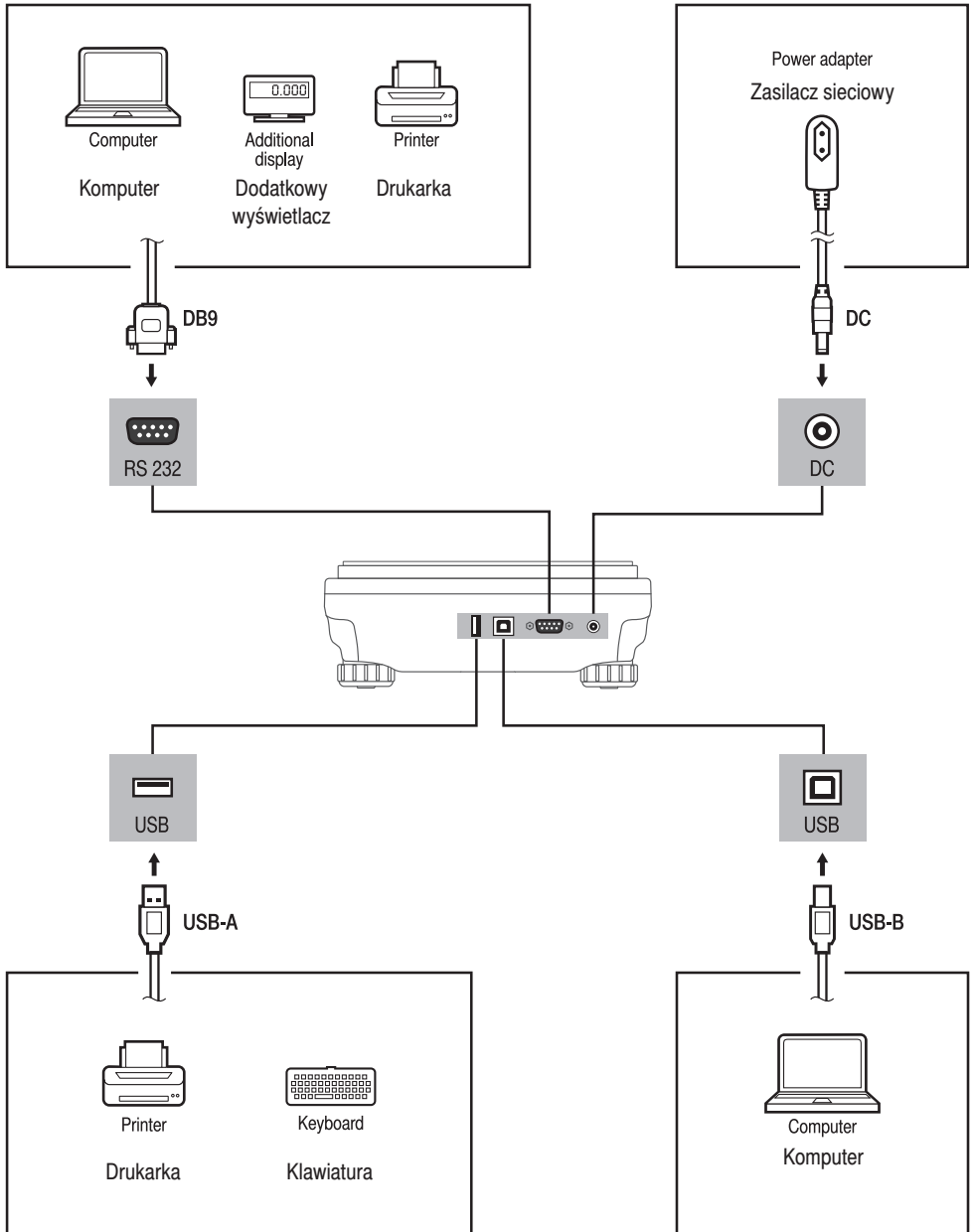
*(lb) pound unit unavailable for verified balances.

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
10. DEVICE CLEANING

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
10.1. General recommendations

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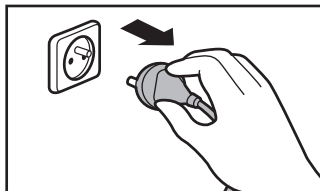
Prior to cleaning it is necessary to disconnect the device from the mains! Remove the plug from the outlet, disconnect all cables.

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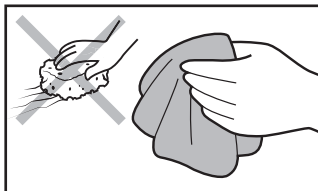Use soft cloths made of microfiber, natural fiber or man-made fiber. Avoid using abrasive cloths or cloths that might scratch the surface.

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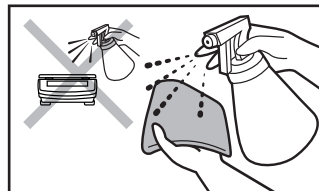Apply the cleanser onto the cloth first. Avoid applying the cleanser directly onto the device.

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Do not use aggressive cleaning agents (e.g. solvents, chlorine preparations, corrosive chemicals, bleach).

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Neither use substances of sharp structure nor intended for scrubbing.

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Make sure that neither dust nor liquid gets inside the weighing system (device inside).

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- **GLASS PANES AND COMPONENTS** can be cleaned using window cleaner.
- **WEIGHING PANS AND STAINLESS STEEL OR ALUMINIUM COMPONENTS** can be cleaned using a cloth slightly soaked in a solution of water and gentle cleanser (e.g. soap or dishwashing liquid).
- **OPERATION PANEL AND HOUSING** can be cleaned using a cloth slightly soaked in a solution of water and gentle cleanser (e.g. soap or dishwashing liquid).
- Dry sample leftovers can be removed using either brush or handheld vacuum cleaner.
- Clean components can be dried using soft cloth or dust-free paper towel, with this the remaining moist will be absorbed.
- **Install all the components providing they are completely dry.**

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